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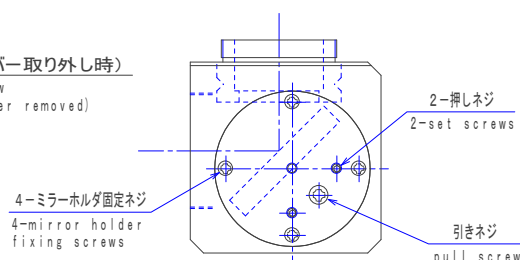
A-A arrow view
(With the cover removed)

I. Overview

By combining with the observation unit with coaxial illumination (OUCI), the observation direction (objective lens mounting direction) can be bent 90 degrees.

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III. Setup

1). Assembling

- Loosen the adapter fixing screws (M3 hexagon socket set screw, 90-degree direction, 2 pieces) and remove the adapter.
- Screw the removed adapter into OUCI, attach the mirror case, and tighten the adapter fixing screw.

2). Adjustment

- Loosen the cover fixing screw and remove the cover.
- You can fine-tune the direction of rotation of the mirror by loosening the mirror holder fixing screws (4 pieces).
- You can fine-tune the tilt (+ direction) of the mirror by loosening the pull screw and tightening the set screw (M2.5 hexagon socket set screw, 2 pieces).
- You can fine-tune the tilt (-direction) of the mirror by loosening the set screw (M2.5 hexagon socket set screw, 2 pieces) and tightening the pull screw.

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IV. Others

- 1). The contents of this manual do not permit the implementation or use of rights such as intellectual property rights owned by SIGMA KOKI Co., Ltd. Reprinting or copying of part or all of this document requires the prior written consent of SIGMA KOKI Co., Ltd.
- 2). The contents of this manual are as of the time of publication and are subject to change without notice.
- 3). The contents of this manual have been carefully prepared, but if you find any mistakes in the description of the materials, please contact the place of purchase or our sales department.
- 4). Combinations other than those described in this manual, or custom specifications may differ from the contents of this manual. If there are products that can be used in combination, also read the instruction manuals for those products.